




Technical Service Consultants Ltd


Microbiology House, Fir St, Heywood, Lancashire OL10 1NW, UK
T : +44 (0)1706 620600 F : +44 (0) 1706 620445
E: sales@tscswabs.co.uk W: www.tscswabs.co.uk

STERILE SAMPLING SPONGE


WITH NEUTRALISING BUFFER

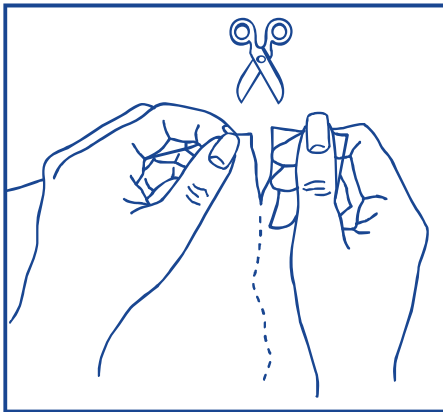


**TEAR, OR CUT OPEN
THE BAG**

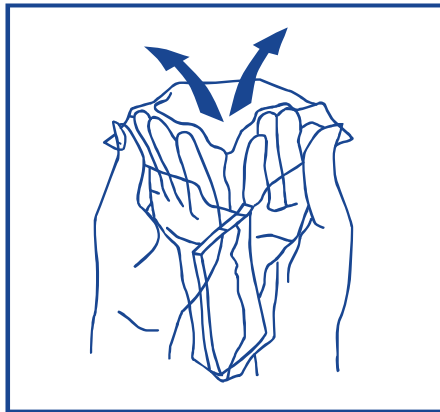


**GRIP EACH SIDE OF THE
BAG WITH FOREFINGER
AND THUMB AND PUSH
THE INSIDE UPWARDS**

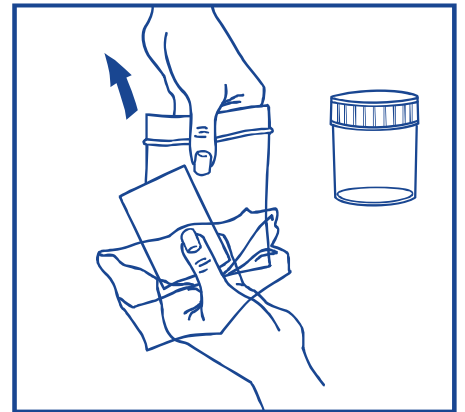


**REMOVE SMALL GRIPPER
BAG, OR TUB, AND PUT TO
ONE SIDE**

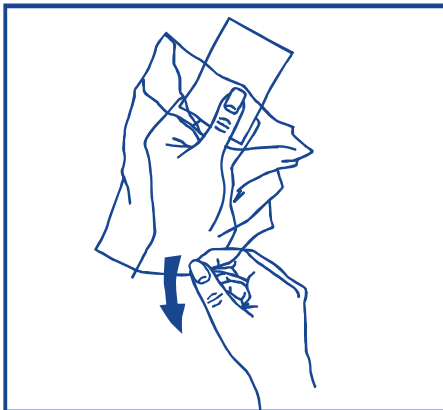

**GRIPPING THE SPONGE
THROUGH THE BAG, PULL
REVERSED BAG OVER HAND**

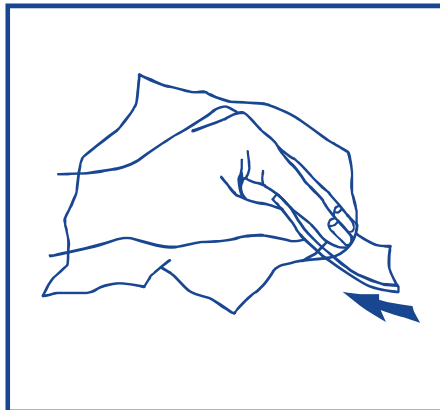

**USE THE SPONGE TO
COLLECT SAMPLE IN
NORMAL WAY**

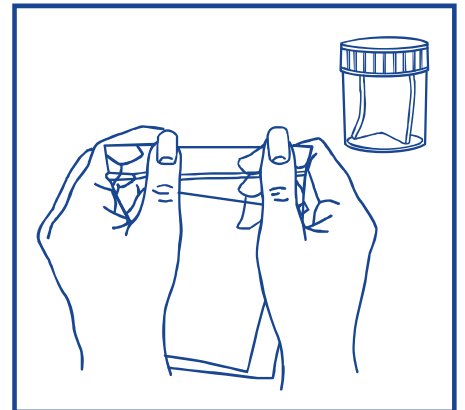

**PLACE THE SPONGE IN THE
GRIPPER BAG, OR TUB,
AND SEAL**

AT NO TIME SHOULD BARE HANDS COME INTO CONTACT WITH THE SPONGE