

GUIDANCE FOR USE


PROTECT Cryopreservation System

All samples should be kept cool until testing is performed.

SETUP FOR FREEZING

STEP 1

Make a suspension of the organism in the vial liquid.

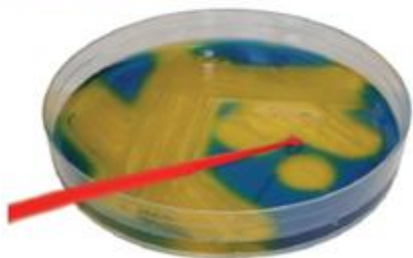

STEP 2

Shake vial to coat the beads.
Stand for 30 seconds.



STEP 3

Remove vial liquid using pastette (TS/126-H).



STEP 4

Label, place in freezer.
Done!



RECOVERY FROM FREEZING

STEP 5

Remove one bead using a sterile needle (TS/3-C).



STEP 6

Roll bead directly onto medium or place in broth and incubate.



Have you considered? Fine Tip Pastettes - TS/126-H



Technical Service
Consultants Ltd

www.tscswabs.co.uk
Tel: +44 (0)1706 620600

GFU_PROTECT_v2 - 30/07/2013